



INSTRUCTION MANUAL



Laboratory Equipment Pty Ltd

Laboratory Equipment Pty Ltd
email: sales@labec.com.au
Ph: 02 9560 2811 • Fax: 02 9560 6131
www.labec.com.au

LABORATORY EQUIPMENT PTY LTD

26 – 30 Farr St, Marrickville, NSW 2204
PHONE: +61 2 95602811 FAX: +61 2 95606131
EMAIL: sales@labec.com.au
www.labec.com.au




INSTRUCTION MANUAL FOR

SI-100

VORTEX MIXER



PLEASE READ THIS MANUAL CAREFULLY BEFORE OPERATION

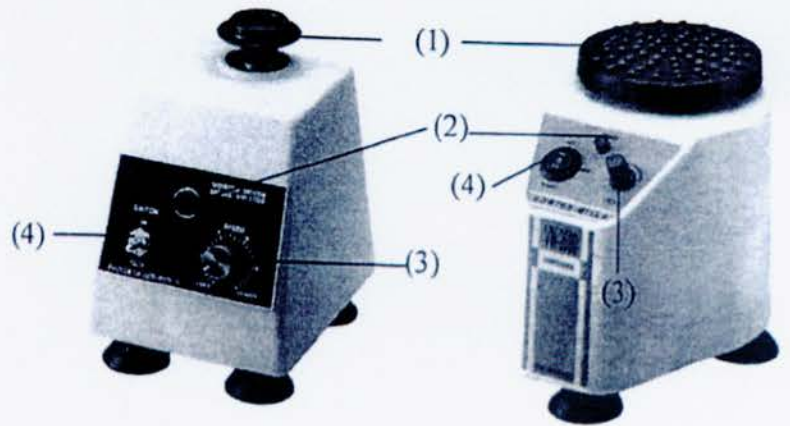

1. RUBBER HEAD
2. POWER LAMP
3. SPEED CONTROL
4. SWITCH

SI-100: ON/ TOUCH

SI-200: AOUTO/OFF/TOUCH

1. Insert the power wire into the socket.
2. Place the "SWITCH" (4) at "TOUCH" Position. The unit will run as soon as touch on the "RUBBER HEAD" (1). The "POWER LAMP" (2) lights.
3. Place the "SWITCH" (4) at the "ON" or ("AUTO" for SI-200) position, the unit will run automatically. The "POWER LAMP" (2) lights.
4. Turn the "SPEED CONTROL" (3) to the required speed.

Warning: Do not make the motor race rotation to avoid damage. After finishing operation, please place the switch at "TOUCH" or "OFF" position.